# Function Check with the ReproEasy Pad

## Quick Reference Guide for Sartorius Models MA37|160


The performance test on models MA37|160 enables you to monitor your device's function, testing both the heating unit and the weighing system at the same time at the temperature you selected.

### 1. Start the performance test


- Select "MA Performance Test" under the "Calibration/Adjustment" item in the Setup menu.
- Select a temperature close to the one you use from the list of temperatures suggested.

### 2. The device now shows you the next work steps

### 3. Place the reference sample on the sample pan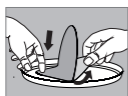

- Open the packaging of the ReproEasy Pad and remove the pad.
- Remove the protective film.

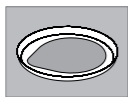- Affix the pad to the center of the sample pan, sticky side down.

### 4. Place the sample pan with the reference sample "ReproEasy Pad" on the pan retainer.

### 5. Close the Hood

- The performance test starts automatically.

### Evaluation of performance test

The device is in a perfect running state and displays "Device is OK" if the Moisture Analyzer performance index (MApi) value falls within the following ranges:

From serial no.: 00343xxxxx

Temperature adjustment

85 °C

125 °C

165 °C

Tolerance range

between 9 and 11 MApi

between 16 and 18 MApi

between 21 and 25 MApi

Up to and including serial no.: 00342xxxxx

Temperature adjustment

70 °C

100 °C

130 °C

Tolerance range

between 9 and 11 MApi

between 16 and 18 MApi

between 21 and 25 MApi

Due to the effect temperature has on the results, the test is only possible once the device has cooled down. If the device is too warm, a relevant message is displayed:

### MA performance test

The device is too warm for a performance test.

The device is still cooling down.

Contact the Sartorius Service Center if the message „Device is not OK.“ appears.

Sartorius Lab Instruments GmbH & Co. KG  
Otto-Brenner-Strasse 20  
37079 Goettingen, Germany  
Phone +49 (0)551 308 0  
www.sartorius.com

Specifications subject to change without notice.  
Copyright Sartorius Lab Instruments GmbH & Co. KG.  
Printed in the EU on paper bleached without chlorine.  
Publication No.: WMA6070-a161102 · Ver. 11 | 2016

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