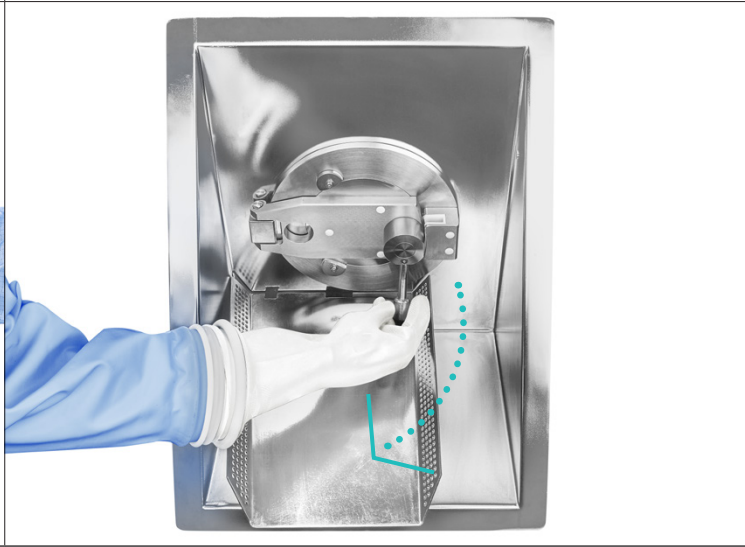


Biosafe® 110 Port

1. Connect the Biosafe® 110 bag	2. Lock	3. Push the lever down to the "Open" position	4. Open the door from the inside, transfer the components, and close the door after the transfer
			
5. Push the lever up to the "Close" position	6. Push the lever down to the "Neutral" position	7. Unlock the Biosafe® 110 bag	8. Remove it
			

Germany


Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen
Phone +49 551 308 0
www.sartorius.com

USA

Sartorius Stedim North America Inc.
565 Johnson Avenue
Bohemia, NY 11716
Toll-Free +1 800 368 7178

Specifications subject to change without notice.
Copyright Sartorius Stedim Biotech GmbH.
DIR No.: 2605342-000-00
Status: 06 | 2020

Biosafe® 110 Port with 90° Lever

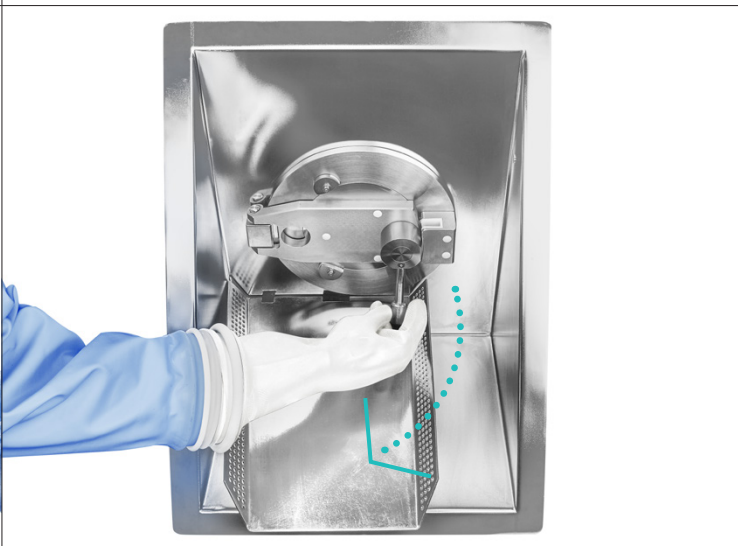1. Connect the Biosafe® 110 bag	2. Lock	3. Push the lever down to the "Open" position	4. Open the door from the inside, transfer the components, and close the door after the transfer
			
5. Push the lever up to the "Close" position	6. Push the lever down to the "Neutral" position	7. Unlock the Biosafe® 110 bag	8. Remove it
			

Germany

Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen
Phone +49 551 308 0
www.sartorius.com

USA

Sartorius Stedim North America Inc.
565 Johnson Avenue
Bohemia, NY 11716
Toll-Free +1 800 368 7178

Specifications subject to change without notice.
Copyright Sartorius Stedim Biotech GmbH.
DIR No.: 2605342-000-00
Status: 06 | 2020